

BAX® System

Quant Solution

BAX® System Quant Solution is a media supplement intended for use with specific quantification protocols. When added to BAX® MP Media, this proprietary solution improves enrichment and recovery, particularly for *Salmonella*.

Features & Benefits:

- Improved post-enrichment recovery of *Salmonella*
- 25 to 50 uses per bottle, depending on protocol
- Liquid format for easier handling compared to other media supplements



How to Use:

BAX System Quant Solution is a clear, colorless, liquid media supplement for specific quantification protocols, particularly for *Salmonella*.

1. Prepare 1 L of BAX MP Media according to the instructions.
2. Cool to 45 to 55 °C.
3. Immediately before use, aseptically transfer 0.5 mL or 1.0 mL of Quant Solution to 1 L of cooled BAX MP Media, and mix.

Note: The volume of Quant Solution added to the BAX MP media depends on your application. For specific protocols, refer to the *BAX System PCR Pathogen Quantification Instructions* (available at www.hygiena.com/documents).

4. Use within 24 hours. Discard remaining media with supplement according to internal SOPs as required by national and local regulations.

Storage & Stability:

Store Quant Solution in its original container at 2 to 8 °C.

Do not use past the expiration date found on the Certificate of Analysis (COA). To find your COA, use the lot number on the bottle and the *COA Search* at www.hygiena.com/documents.

Product No.	Description	Quantity
MED2032	BAX® System Quant Solution	25 mL
MED2003	BAX® System MP Media	2.5 kg
MED2029	BAX® System MP Media	10 kg